## USB DIGITAL MICROSCOPE

GB

### USB DIGITAL MICROSCOPE

Operation and safety notes

DE AT CH

### USB- DIGITALMIKROSKOP

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**A**

<b>Warnings and symbols used</b> .....	Page	6
<b>Introduction</b> .....	Page	7
Intended use .....	Page	7
Trademark notices .....	Page	7
Scope of delivery .....	Page	7
Parts list .....	Page	7
Technical data .....	Page	8
<b>Safety instructions</b> .....	Page	9
<b>Before first use</b> .....	Page	11
Unpacking the product .....	Page	11
Downloading the software .....	Page	12
Setting up the product .....	Page	12
<b>Use</b> .....	Page	12
Adjusting the brightness of the LED ring light .....	Page	13
Adjusting the stand .....	Page	13
Adjusting the zoom .....	Page	13
Taking a photo .....	Page	13
Software overview for Windows® computers .....	Page	13
Software overview for MacOS computers .....	Page	17
<b>Troubleshooting</b> .....	Page	20
<b>Cleaning and care</b> .....	Page	21
Cleaning .....	Page	21
Storage .....	Page	21
<b>Disposal</b> .....	Page	22
<b>Warranty</b> .....	Page	22
Warranty claim procedure .....	Page	23
<b>Service</b> .....	Page	24

## Warnings and symbols used

The following warnings are used in this user manual and on the packaging:

	<p><b>WARNING!</b> This symbol in combination with the signal word "Warning" marks a medium-risk hazard that if not prevented could result in death or serious injury.</p>		Warning against optical radiation
			Use the product in dry indoor spaces only.
	<p><b>CAUTION!</b> This symbol in combination with the signal word "Caution" marks a low-risk hazard that if not prevented could result in minor or moderate injury.</p>		Direct current/voltage
			LED button <span style="border: 1px solid black; padding: 0 2px;">8</span>
			Photo button <span style="border: 1px solid black; padding: 0 2px;">7</span>
	<p><b>ATTENTION!</b> This symbol with the signal word "Attention" indicates a possible property damage.</p>		Safety information Instructions for use
	<p><b>NOTE:</b> This symbol in combination with "Note" provides additional useful information.</p>		CE mark indicates conformity with relevant EU directives applicable for this product.

# **USB DIGITAL MICROSCOPE**

## **● Introduction**

We congratulate you on the purchase of your new product. You have chosen a high quality product. The instructions for use are part of the product. They contain important information concerning safety, use and disposal. Before using the product, please familiarise yourself with all of the safety information and instructions for use. Only use the product as described and for the specified applications. If you pass the product on to anyone else, please ensure that you also pass on all the documentation with it.

## **● Intended use**

- This product is used to magnify small objects, and to display images on a computer screen.
- Any other usage or modification of the product is deemed to be improper and carries a significant risk of accidents.
- The manufacturer accepts no responsibility for damages resulting from improper usage.

- This product is not intended for commercial use.
- Use the product in dry indoor spaces only.

## **● Trademark notices**

- USB® is a registered trademark of USB Implementers Forum, Inc.
- Windows® is a registered trademark of Microsoft Corporation in the United States and other countries.
- Any other names and products are trademarks or registered trademarks of their respective owners.

## **● Scope of delivery**

- 1 USB digital microscope
- 1 Stand
- 1 Storage case
- 1 User manual

## **● Parts list**

Fig. A

- 1 Clamp
- 2 Bracket
- 3 Locking screw
- 4 Stand
- 5 Storage case

- 6 Digital microscope
- 7  button (Photo button)
- 8  button (LED button)
- 9 Zoom wheel
- 10 LED ring lighting
- 11 Protection cap
- 12 Objective lens
- 13 USB cable
- 14 USB plug (Type C)
- 15 USB adapter (Type A)

## ● Technical data

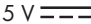
Required connections:	1 USB port
Operating voltage:	5 V 
Current consumption:	216 mA
Cable length:	approx. 1.36 m
Operating systems:	Windows® 8 or above macOS X V10.5 or above
Video format:	Windows®: AVI (up to 30 fps) MacOS: MP4 (up to 30 fps)
Image format:	PNG, JPG

Image resolution:	640 × 480 pixels/ 1,280 × 720 pixels */ 1,920 × 1,080 pixels *
Magnification:	50× to 1,000×
Working distance:	15 mm (1,000×) to 40 mm (50×)
Illumination:	8 LEDs
Computer connection:	USB Type A ** USB Type C **
Operating temperature:	+5 °C to +35 °C
Storage temperature:	0 °C to +45 °C
Storage humidity:	10 % to 70 % relative humidity
Altitude:	Max. 2,000 m above sea level
Pollution degree :	2
Dimensions (L × W × H):	approx. Ø 32 × 120 mm
Weight:	approx. 171 g
* Interpolated	
** Data transfer and power supply are carried out via the USB port on the computer at the same time.	

- The specifications and design are subject to change without prior notification.

## Safety instructions

**BEFORE USING THE PRODUCT, FAMILIARISE YOURSELF WITH ALL OF THE SAFETY INFORMATION AND INSTRUCTIONS FOR USE! WHEN PASSING THIS PRODUCT ON TO OTHERS, ALSO INCLUDE ALL THE DOCUMENTS!**

In the case of damage resulting from non-compliance with these operating instructions the warranty claim becomes invalid!

No liability is accepted for consequential damage! In the case of material damage or personal injury caused by incorrect handling or non-compliance with the safety instructions, no liability is accepted!

**⚠ WARNING! DANGER TO LIFE AND RISK OF ACCIDENT FOR INFANTS AND CHILDREN!**

- Never leave children unsupervised with the packaging material. The packaging material represents a danger of suffocation.

Children frequently underestimate the dangers. Always keep children away from the packaging material.

- This product contains small parts that can be swallowed by children. If small parts are swallowed, consult a doctor immediately.
- This product may be used by children aged 8 years and above, as well as by persons with reduced physical, sensory or mental capacities or lacking experience and/or knowledge, as long as they are supervised or instructed in the safe use of the product and understand the associated risks. Do not allow children to play with the product. Cleaning and user maintenance should not be performed by children without supervision. The product is not a toy.

- ⚠ WARNING! Risk of injury!** This product contains components and/ or accessories that can cause minor or serious body injury if

used improperly. It is therefore essential that you observe the following safety information in order to prevent body injuries.

- Sharp-edged and pointed tools are often used for working with this product. Keep this product and all accessories and tools in a place inaccessible to children.
- Chemicals and liquids do not belong in children's hands! Do not drink! Clean hands thoroughly under running water after use. In case of accidental contact with eyes or mouth, rinse with water. In case of complaints, consult a doctor immediately and show the substances.
- Inspect the product before each use! Discontinue use if there are any signs of damage.
- Sudden temperature changes may cause condensation to form inside the product. To prevent short circuits, allow the product to acclimatise to its environment before using it again.
- Do not use the product near heat sources such as radiators or other heat-emitting devices!
- Do not throw the product into fire or expose it to high temperatures.
- Never open the product! It contains no internal parts that require maintenance.
- Do not expose the product to direct heat sources (e.g. heating).
- Avoid any contact with splashing or dripping water and corrosive liquids, and do not use the product in the vicinity of water. In particular, the product should never be submerged in liquid. Do not place any objects filled with liquid (e.g. vases or drinks) on or near the product.
- Do not use the product in the immediate vicinity of magnetic fields (e.g. speakers).
- Do not place sources of fire (e.g. burning candles) on or near the product.
- Do not insert any objects into the interior of the product.

- Do not position the equipment so that it is difficult to operate the USB port.
- Do not subject the product to excessive shocks and vibrations.
-  Do not look directly into the light beam or look at the light beam with optical devices.
- OWIM GmbH & Co KG shall not be liable for interference with radios or televisions due to unauthorised modification of the product.
- The user of the product is fully responsible for correcting interference caused by such unauthorised modification of the product, as well as replacement of such products.
- The product must be repaired if it has been damaged in any way (for example, if the housing is damaged, if liquid or objects have penetrated the product or if the product has been exposed to rain or moisture).
- A repair is also required if the product is not working properly or has been dropped. If smoke is produced, or there are any unusual sounds or smells, disconnect the product from the computer. In such cases, the product must not be used until it has been inspected by authorised service personnel. The product must only be repaired by qualified personnel. Never open the housing of the product.

## ● **Before first use**

### ● **Unpacking the product**

1. Take the product out of the packaging and remove all packaging materials.
2. Check to make sure that all listed parts are included (see “Scope of delivery”).
3. Check whether the product and all parts are in good condition, if any damage or defect is detected, do not use the product, but follow the procedure described in chapter “Warranty”.

## ● Downloading the software

- ① **NOTE:** Use the **HiView** software to view, save and label the images of the digital microscope **6**.

### For Windows® computers

1. Open the website [www.lidl-service.com](http://www.lidl-service.com).  
The QR code is shown on the back page of the user manual.
2. Searching for the software:  
Enter the product number IAN 490136\_2501.
3. Download the **HiView** software to your computer.

### For MacOS computers

1. Open the App Store and search for **HiView**.
2. Download the **HiView** software to your computer.

## ● Setting up the product

1. Snap the digital microscope **6** into the clamp **1**.
2. Switch on the computer.

3. Optional: Select your preferred USB plug type by attaching or detaching the USB adapter (Type A **15**) on the USB plug (Type C **14**).
4. Connect the USB plug (Type C **14** or Type A **15**)
  - to a free USB port on your computer
  - or
  - to a USB hub that is connected to your computer.The LED ring lighting **10** lights up when the USB plug (Type C **14** or Type A **15**) successfully connect to the computer.

5. Remove the protection cap **11** from the digital microscope **6**.
6. Place the observation object at the working distance (see “Technical data”) below the objective lens **12**.

## ● Use

- ① **NOTE:** If the image does not appear sharp on the computer screen: Carry out the following steps again:
- Adjusting the brightness of the LED ring light

- Adjusting the stand
- Adjusting the zoom

## ● Adjusting the brightness of the LED ring light

- Press the  button **8** once or multiple times, until the LED ring light **10** illuminates the observation object well.

## ● Adjusting the stand

1. Loosen the locking screw **3**.
2. Select the required angle of the stand.
3. Tighten the locking screw **3**.

## ● Adjusting the zoom

- Slowly rotate the zoom wheel **9** towards 50× or 1,000×, until the observation object is clearly visible. The image appears sharp on the computer screen.

## ● Taking a photo

1. To avoid motion blur: Keep the digital microscope **6** steady.
2. Press the  button **7**.  
or  
Tap  on the graphic user interface of **HiView**.



## ● Software overview for Windows® computers

### Main interface

The following settings are available in the main interface:


Interface	Function
Name: <input type="text" value="USB 2.0 Camera"/> 	<input type="checkbox"/> Select your digital microscope
Resolution: <input type="text" value="640x480"/> 	<input type="checkbox"/> Set the resolution
 Photo	<input type="checkbox"/> Take a photo

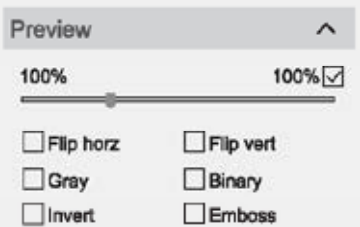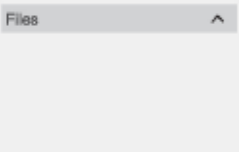Interface	Function
 <b>Video</b>	<ul style="list-style-type: none"> <li><input type="checkbox"/> Record a video</li> </ul>
 <b>Files</b>	<ul style="list-style-type: none"> <li><input type="checkbox"/> Open the folder that contains the recorded images</li> <li><input type="checkbox"/> The designated folder can be changed under  <b>Setting</b></li> </ul>
 <b>Setting</b>	<ul style="list-style-type: none"> <li><input type="checkbox"/> Select the appearance: <b>Ruler/Line/Text</b></li> <li><input type="checkbox"/> Select the folder in which the recorded images are saved</li> <li><input type="checkbox"/> Select the file format for images</li> <li><input type="checkbox"/> Change the on-screen language</li> <li><input type="checkbox"/> For further details: see "Setting"</li> </ul>

Interface	Function
	<ul style="list-style-type: none"> <li><input type="checkbox"/> Zoom in/out on the image</li> <li><input type="checkbox"/> Picture effects: <ul style="list-style-type: none"> <li>– <b>Flip horz:</b> Flip the image horizontally</li> <li>– <b>Flip vert:</b> Flip the image vertically</li> <li>– <b>Grey:</b> Convert colour into grey scale</li> <li>– <b>Binary:</b> Convert colour into black and white</li> <li>– <b>Invert:</b> Invert the colour on the image (Artistic effect)</li> <li>– <b>Emboss:</b> Embossing (Artistic effect)</li> </ul> </li> </ul>
	<ul style="list-style-type: none"> <li><input type="checkbox"/> Recorded images and videos</li> </ul>

## Setting Setting

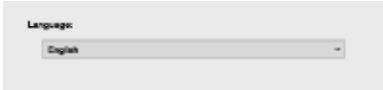The following settings are available:

Interface	Function
	<ul style="list-style-type: none"> <li><input type="checkbox"/> Select the colour of the ruler</li> <li><input type="checkbox"/> Select the colour and size of the line and text on the image</li> </ul>
	<ul style="list-style-type: none"> <li><input type="checkbox"/> Select the folder for saving the photos and videos</li> <li><input type="checkbox"/> Select the format for saving the image</li> </ul>
	<ul style="list-style-type: none"> <li><input type="checkbox"/> Select the language of the software</li> <li><input type="checkbox"/> Available languages: English, German, French, Spanish, Italian, Portuguese</li> </ul>

## Labelling tools

The following tools are available:

Tool	Function	Tool	Function
	<ul style="list-style-type: none"> <li><input type="checkbox"/> Showing the cross hairs for reference</li> </ul>		<ul style="list-style-type: none"> <li><input type="checkbox"/> Draw and measure an angle</li> </ul>

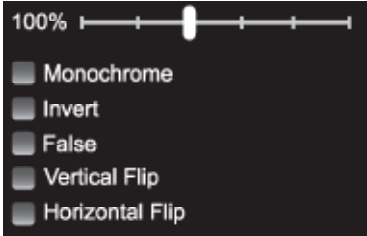
Tool	Function	Tool	Function
	<input type="checkbox"/> Draw a line		<input type="checkbox"/> Arbitrary draw
	<input type="checkbox"/> Draw a parallel line		<input type="checkbox"/> Add text
	<input type="checkbox"/> Draw a rectangle		<input type="checkbox"/> Move the image
	<input type="checkbox"/> Draw a circle		<input type="checkbox"/> Return to previous operation
	<input type="checkbox"/> Draw a concentric circle		<input type="checkbox"/> Delete all annotations

## ● Software overview for MacOS computers

### Main interface

The following settings are available in the main interface:

Interface	Function
	<input type="checkbox"/> Select your digital microscope 
	<input type="checkbox"/> Set the resolution
	<input type="checkbox"/> Take a photo
	<input type="checkbox"/> Record a video

Interface	Function
	<ul style="list-style-type: none"> <li><input type="checkbox"/> Open the folder that contains the recorded images</li> <li><input type="checkbox"/> The designated folder can be changed under </li> </ul>
	<ul style="list-style-type: none"> <li><input type="checkbox"/> Select the appearance: <b>Line/Text</b></li> <li><input type="checkbox"/> Select the folder in which the recorded images are saved</li> <li><input type="checkbox"/> Select the file format for images</li> <li><input type="checkbox"/> For further details: see "Setting"</li> </ul>
	<ul style="list-style-type: none"> <li><input type="checkbox"/> Zoom in/out on the image</li> <li><input type="checkbox"/> Picture effects: <ul style="list-style-type: none"> <li>– <b>Monochrome:</b> Turn colour into grey scale</li> <li>– <b>Invert:</b> Invert the colour on the image (artistic effect)</li> <li>– <b>False:</b> False colour (artistic effect)</li> <li>– <b>Vertical Flip:</b> Flip the image vertically</li> <li>– <b>Horizontal Flip:</b> Flip the image horizontally</li> </ul> </li> </ul>

Interface	Function
	<input type="checkbox"/> Recorded images and videos

## Setting

The following settings are available:

Interface	Function
	<input type="checkbox"/> Select the colour and size of the line and text on the image
	<input type="checkbox"/> Select the folder for saving the photos and videos <input type="checkbox"/> Select the format for saving the image

## Labelling tools

The following tools are available:

Tool	Function	Tool	Function
	<input type="checkbox"/> Showing the cross hairs for reference		<input type="checkbox"/> Arbitrary draw

Tool	Function	Tool	Function
	<input type="checkbox"/> Draw a line		<input type="checkbox"/> Move the image
	<input type="checkbox"/> Draw a parallel line		<input type="checkbox"/> Add an arrow
	<input type="checkbox"/> Draw a circle		<input type="checkbox"/> Add text
	<input type="checkbox"/> Draw a concentric circle		<input type="checkbox"/> Return to previous operation
	<input type="checkbox"/> Draw a rectangle		<input type="checkbox"/> Delete all annotations
	<input type="checkbox"/> Draw and measure an angle		

## ● Troubleshooting

Problem	Possible cause	Solution
The product does not function.	The USB plug <a href="#">14</a> or the USB adapter <a href="#">15</a> are not properly connected to the computer.	Remove the USB plug <a href="#">14</a> or the USB adapter <a href="#">15</a> and reconnect to the computer again.
	The protection cap <a href="#">11</a> covers the digital microscope <a href="#">6</a> .	Remove the protection cap <a href="#">11</a> from the digital microscope <a href="#">6</a> .

Problem	Possible cause	Solution
The image does not appear sharp on the computer screen.	The zoom is not adjusted properly.	Adjust the zoom (see "Adjusting the Zoom").
	The observation object is too close or too far away from the digital microscope <b>6</b> .	Adjust the position of the observation object.
	There is motion blur on the images.	When pressing the  button <b>7</b> : Keep the digital microscope <b>6</b> steady.

## ● Cleaning and care

### ● Cleaning

#### ① **NOTE:**

- Do not use chemical, alkaline, abrasive or other aggressive detergents or disinfectants to clean this product as they might be harmful to its surfaces.
- Never allow fluids to get into the product.
- Before cleaning: Attach the protection cap **11** to the digital microscope **6**.

- Clean the product with a soft, dry cloth.

### ● Storage

- Clean the product (see "Cleaning").
- Insert the product and its accessories into the storage case **5**. Store the storage case in a dry indoor area out of direct sunlight.
- Store the product in a place that is inaccessible to children.

## ● Disposal

### **Packaging:**

The packaging is made of environmentally friendly materials, which you be disposed through your local recycling facilities.

Observe the marking of the packaging materials for waste separation, which are marked with abbreviations (a) and numbers (b) with following meaning: 1-7: plastics/20-22: paper and fibreboard/80-98: composite materials.

### **Product:**

Contact your local refuse disposal authority for more details of how to dispose of your worn-out product.

The adjacent symbol of a crossed out dustbin on wheels indicates this product is subject to directive 2012/19/EU. This directive states at the end of the life this product

must not be disposed of through regular household refuse but must be returned to special collection sites, recycling depots or waste management companies.

### **This disposal is free of charge to you.**

Protect the environment and dispose of properly.

## ● Warranty

The product has been manufactured to strict quality guidelines and meticulously examined before delivery. In the event of material or manufacturing defects you have legal rights against the retailer of this product. Your legal rights are not limited in any way by our warranty detailed below.

The warranty for this product is 5 years from the date of purchase. The warranty period begins on the date of purchase. Keep the original sales receipt in a safe location as this document is required as proof of purchase.

Any damage or defects already present at the time of purchase must be reported without delay after unpacking the product.

Should the product show any fault in materials or manufacture within 5 years from the date of purchase, we will repair or replace it – at our choice – free of charge to you. The warranty period is not extended as a result of a claim being granted. This also applies to replaced and repaired parts.

This warranty becomes void if the product has been damaged, or used or maintained improperly.

The warranty covers material or manufacturing defects. This warranty does not cover product parts subject to normal wear and tear, thus considered consumables (e.g. batteries, rechargeable batteries, tubes, cartridges), nor damage to fragile parts, e.g. switches or glass parts.

## ● **Warranty claim procedure**

To ensure quick processing of your claim, observe the following instructions:

Make sure to have the original sales receipt and the item number (IAN 490136\_2501) available as proof of purchase.

You can find the item number on the rating plate, an engraving on the product, on the front page of the instruction manual (bottom left), or as a sticker on the rear or bottom of the product.

If functional or other defects occur, contact the service department listed below either by telephone or by e-mail.

Once the product has been recorded as defective you can return it free of charge to the service address that will be provided to you. Ensure to enclose the proof of purchase (sales receipt) and a short, written description outlining the details of the defect and when it occurred.

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**OWIM GmbH & Co. KG**